

STANDARD OPERATING PROCEDURES

DIVISION OF COMPARATIVE MEDICINE
UNIVERSITY OF SOUTH FLORIDA

SOP#: 1146

Date Issued: 3/07

Date Revised: New

Page 1 of 2

TITLE:	PRL Power Foamer
SCOPE:	Animal Care Personnel
RESPONSIBILITY:	Facility Manager and Technical Staff
PURPOSE:	To Outline the Proper Procedures for Use and Maintenance of a PRL Power Foamer

I. PURPOSE

1. This procedure outlines the use and maintenance of the PRL Power Foamer used for the application of a chemical solution (i.e., Virkon[®]S) in the form of thick foam.

II. RESPONSIBILITY

1. It is the responsibility of the Facility Manager to ensure that portable equipment is appropriately maintained in good working order.
2. It is the responsibility of the Facility Manager to ensure that technical staff using this equipment are adequately trained and experienced in the use of the PRL Power Foamer.
3. It is the responsibility of the technical staff using this equipment to read the equipment manual and understand the operation of the PRL Power Foamer and associated chemical use.

III. EQUIPMENT USE

1. Foaming, when done properly, reduces the likelihood of chemical splashes, the creation

