

# STANDARD OPERATING PROCEDURES

DIVISION OF COMPARATIVE MEDICINE  
UNIVERSITY OF SOUTH FLORIDA

SOP#: 1161.2

Date Issued: 5/15

Date Revised: 7/20

Page 1 of 2

---

TITLE:	Germinator™ 500 Glass Bead Dry Sterilizer
SCOPE:	Animal Care Personnel
RESPONSIBILITY:	Facility Manager, Professional and Administrative Staff
PURPOSE:	To Outline the Proper Procedures for Use and Maintenance of a Germinator™ Glass Bead Sterilizer

---

## I. PURPOSE

1. This procedure outlines the use and maintenance of the Germinator™ Glass Bead Sterilizer used to

